

## CellPlayer™ NuLight Mix: HT-1080 NuLight Green and A549 NuLight Red

EsSEN BioScience Catalog Number: 4516

### Storage

Liquid Nitrogen

*Note: Cells can be thawed and cultured immediately upon receipt or stored in liquid nitrogen for long-term storage. Storage at -80°C is not recommended. Cells should be used within 1 year of delivery.*

### Presentation

1mL,  $1 \times 10^6$  cells/mL in 90% FBS, 10% DMSO

- HT-1080 NuLight Green:  $5 \times 10^5$  cells/vial
- A549 NuLight Red:  $5 \times 10^5$  cells/vial

### Recommended Media and Components

Ham's F-12 + GlutaMAX™-1 (Cat# 31765 Gibco/Life Technologies; Low riboflavin media for fluorescence imaging)



10% FBS (Cat# SH30071 Thermo Hyclone)

1% Pen/Strep (Cat# 15140 Gibco/Life Technologies)

0.5µg/mL Puromycin (Cat# A11138-03 Gibco/Life Technologies)

### Background

Each vial contains a mixed stable population of 500,000 HT-1080 NuLight Green cells (Cat# 4486) and 500,000 A549 NuLight Red cells (Cat# 4491). Parental HT-1080 and A549 cells were purchased from ATCC (Cat# CCL-121 and Cat# CCL-185, respectively). HT-1080 and A549 cells were made individually using the Essen CellPlayer NuLight Green Lentivirus (Cat# 4475; EF1α, puromycin) and NuLight Red Lentivirus (Cat# 4476; EF1α, puromycin), respectively, at an MOI of 3 (TU/cell) in the presence of 8µg/mL polybrene following the standard Essen protocol. This resulted in  $\geq 70\%$  transduction efficiency. Individual cell lines were mixed 1:1 to a total final concentration of 1 million cells/mL in 90% FBS and 10% DMSO. All individual cell lines in our catalog have been certified mycoplasma free by ATCC and our stable populations have been unambiguously authenticated using ATCC's Short Tandem Repeat (STR) profiling.

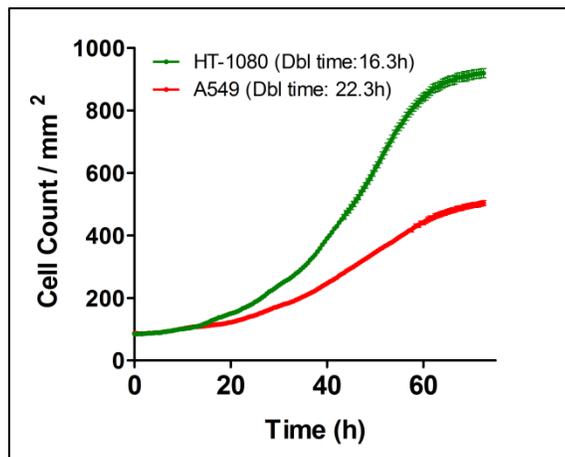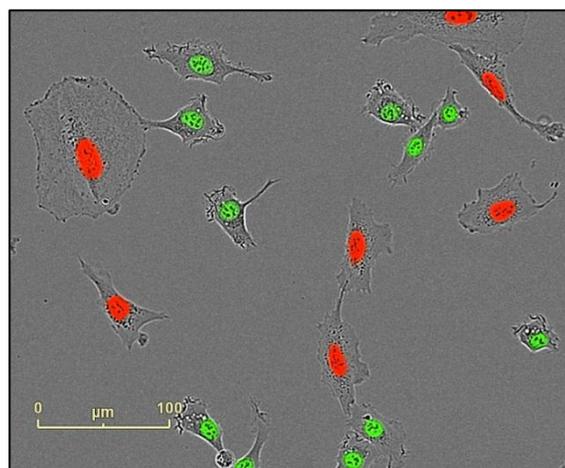

**Figure 1.** (Top) HD-Phase and fluorescence blend of HT-1080 NuLight Green and A549 NuLight Red cells imaged at 20x in IncuCyte ZOOM. (Bottom) Proliferation and doubling times of each cell type in co-culture (4x).



## Protocols and Procedures

### Thawing and Culturing Cells

1. Immediately thaw vial of NuLight Mix cells, or store in liquid nitrogen upon receipt.
2. In a 37°C water bath, quickly thaw each vial by gentle agitation. Be careful not to submerge entire vial to avoid contamination. This process should take no more than 2 minutes. Remove vial when only a tiny ice crystal remains.
3. Wipe vial with 70% ethanol and perform step 4 or 5 in a biosafety hood.
4. Prepare a T75 Flask
  - a. Dispense 5mL culture media into a T25 flask.
  - b. Aliquot cells into the T25 flask.
  - c. Gently rock the culture flasks to evenly distribute the cells and place in 37°C incubator.
  - d. 5-8hr later, replace media containing DMSO with fresh media.
5. Prepare a Microplate
  - a. Dispense 5-10mL media into a 15mL conical tube
  - b. Aliquot cells into the 15mL conical tube and centrifuge at 1,000rpm for 5 minutes.
  - c. Aspirate media, re-suspend the cell pellet in fresh media, and distribute to wells according to Table 1.
  - d. Let cells sit at room temperature for a minimum time of 30 minutes before incubating at 37°C.

**Table 1: Recommended volumes and final cell densities for microplates**

Microplate Vessel	Volume to re-suspend cell pellet	Volume to add per well	Final cell density per well
6-well	15mL	2.3mL	~150,000 cells/well
12-well	13mL	1mL	~75,000 cells/well
24-well	13mL	0.5mL	~35,000 cells/well
48-well	13mL	0.25mL	~20,000 cells/well
96-well	12mL	0.1mL	~8,000 cells/well

### IncuCyte ZOOM™ Set-up and Data Analysis

1. Immediately place the flask or microplate in IncuCyte ZOOM™. Let the flask or plate warm up for 15-30 minutes.
2. Set the Scan Schedule:
  - a. Channel Selection: Phase, Green, Red
  - b. Spectral Unmixing: 8% of Red removed from Green
  - c. Set interval every 2 hours
  - d. Job Type: Basic Analyzer
  - e. Processing Definition: Phase/Green/Red Processor (4x), (10x), or (20x)
3. Scan for 3-5 days, collect 4-6 images representative of the mixed culture
  - a. Create a processing definition and launch an analysis job following the Fluorescence Processing Technical Note.
  - b. Analyze data following the Proliferation Application Note.





## Related Products

### **NuLight/CytoLight Reagents:**

Cat.# 4475 CellPlayer NuLight Green (Lenti, EF-1 alpha, puro)  
 Cat.# 4481 CellPlayer CytoLight Green (Lenti, EF-1 alpha, puro)  
 Cat.# 4513 CellPlayer CytoLight Green (Lenti, CMV, no selection)

Cat.# 4476 CellPlayer NuLight Red (Lenti, EF-1 alpha, puro)  
 Cat.# 4482 CellPlayer CytoLight Red (Lenti, EF-1 alpha, puro)

### **NuLight Cell Lines:**

Cat.# 4485 CellPlayer HT-1080 NuLight Red  
 Cat.# 4487 CellPlayer MDA-MB-231 NuLight Red  
 Cat.# 4489 CellPlayer HeLa NuLight Red  
 Cat.# 4491 CellPlayer A549 NuLight Red  
 Cat.# 4506 CellPlayer HUVEC NuLight Green  
 Cat.# 4453 CellPlayer HUVEC CytoLight Green

Cat.# 4486 CellPlayer HT-1080 NuLight Green  
 Cat.# 4488 CellPlayer MDA-MB-231 NuLight Green  
 Cat.# 4490 CellPlayer HeLa NuLight Green  
 Cat.# 4492 CellPlayer A549 NuLight Green  
 Cat.# 4511 CellPlayer Neuro-2a NuLight Green  
 Cat.# 4512 CellPlayer Neuro-2a NuLight Red

For additional information on this and other products, please contact Essen BioScience at: [sales@essenbio.com](mailto:sales@essenbio.com).

**For research use only. Not for therapeutic or diagnostic use.**

## Licenses and Warranty

Essen BioScience warrants that this product performs as described on the product label and in all literature associated with the sale of said product when used in accordance with the described protocol. This limited warranty is the sole warranty. No other warranties exist that extend beyond this warranty, either expressed or implied. Essen BioScience disclaims any implied warranty of merchantability or warranty of fitness for a particular purpose. Essen BioScience disclaims any responsibility for injury or damage and shall not be liable for any proximate, incidental or consequential damages in connection with this product.

If it is proven to the satisfaction of Essen BioScience that this product fails to meet performance specifications, Essen BioScience's sole obligation, at the option of Essen BioScience, is to replace the product or provide the purchaser with credit at or below the original purchase price. This limited warranty does not extend to anyone other than the purchaser. Notice of suboptimal performance must be made to Essen BioScience within 30 days of receipt of the product.

This Essen BioScience product contain proprietary nucleic acid(s) coding for proprietary fluorescent protein(s) being, including its derivatives or modifications, the subject of pending patent applications and/or patents owned by Evrogen JSC (hereinafter "Evrogen Fluorescent Proteins").

The purchase of Essen BioScience products incorporating these fluorescent proteins conveys to the buyer the non-transferable right to use Evrogen Fluorescent Proteins only for research conducted by the buyer. No rights are conveyed to modify or clone the gene encoding fluorescent protein contained in this product or to use Evrogen Fluorescent Proteins for commercial purposes. The right to use Evrogen Fluorescent Proteins specifically excludes the right to validate or screen compounds for commercial purposes. For information on commercial licensing, contact Evrogen Licensing Department, email: [license@evrogen.com](mailto:license@evrogen.com).

